



**Instruction Manual
for
Filtration Centrifuge
Sieva-3**

1. GENERAL INFORMATION.....	1
1.1 Danger, precautions and warranty.....	1
1.2 Description.....	2
1.3 Delivery Package.....	2
1.4 Warranty.....	2
1.5 Operating and Display Elements	3
1.5.1 LED light.....	3
1.5.2 LCD-Display	4
1.6 Signs and Indications of the Centrifuge	4
1.6.1 General.....	4
1.6.2 Product Nameplate (Example)	5
1.6.3 Warning and Information Signs	6
2. INSTALLATION	7
2.1 Unpackaging.....	7
2.2 Space Requirements	7
2.3 Installation.....	7
2.4 Installation of the rotor	8
2.4.1 Mounting of the centrifugal basket.....	8
2.4.2 Überladung von Rotoren	8
2.4.3 Removing of the rotor	8
2.4.4 Functional principle of the baskets	9
3. BASIC ADJUSTMENTS	10
3.1 Adjustment of the Rotor Type	10
3.2 Access to the Mode: "Standard settings"	10
3.3 Acoustic Signal Turn On / Off	11
3.4 Volume Pre-Selection of Sound Signal.....	11
3.5 Song Selection – End of Run.....	12
3.6 Keyboard Sound Turn On / Off	12
3.7 Sleep Mode Setting	13
3.8 Retrieving Operation Data	13
4. OPERATION	14
4.1 Power Switch.....	14
4.2 Centrifuge Lid	14
4.2.1 Lid Release	14
4.2.2 Lid Lock.....	15
4.3 Pre-Selection	15
4.3.1 Pre-Selection of Speed and RCF-Value.....	15
4.3.2 Pre-Selection of Running Time.....	16
4.3.3 Pre-Selection of Brake Intensity and Acceleration.....	16
4.4 Program.....	17

4.4.1	Storage of Programs	17
4.4.2	Recall of Stored Programs	18
4.4.3	Leaving Program Mode	19
4.5	Starting and Stopping the Centrifuge	19
4.5.1	Starting the Centrifuge	19
4.5.2	Stopping the Centrifuge	19
4.6	Imbalance Detection	20
4.7	Volume Flow	20
4.8	External Temperature Control	20
5.	MAINTENANCE	21
5.1	Maintenance and Cleaning	21
	Cleaning and Disinfecting of the Unit	22
	Cleaning and Disinfection of the Rotor	22
	Disinfection of Aluminum Rotors	22
	Disinfection of PP-Rotors	23
	Glass Breakage	23
	Lifetime of Rotors, Round and Rectangular Buckets, Accessories	23
6.	TROUBLE SHOOTING	24
6.1	Error Messages: Problem / Solution	24
6.2	Survey of Possible Error Messages and Solutions	24
6.2.1	Lid Release during Power Failure (Emergency Lid Release)	24
6.2.2	Description of the Error Message System	25
6.2.3	Procedure while error 14	25
7.	TRANSPORT, STORAGE AND DISPOSAL	26
	Transport	26
	Storage	26
	Disposal	26
8.	APPENDIX	I
	EC – Conformity Declaration	II
	Tabelle 1: Technical Data	III
	Table 2: Max. Speed and RCF-Values for Permissible Rotors	IV
	Table 3: Permissible Net Weight	IV
	Table 4: Error Messages	V
	Table 5: Abbreviations	V
	Redemption Form / Decontamination Certificate	1

1. GENERAL INFORMATION

1.1 Danger, precautions and warranty

Before putting the centrifuge into operation, please read this instruction manual carefully.



This device may only be operated by trained specialist staff. They must have carefully read the operating manual and be familiar with the function of the device.

To protect people and environment the following precautions should be observed:

- During centrifugation, the presence of people and the setting up of hazardous materials is prohibited within 30 cm around the centrifuge according to the regulations of EN 61010-2-020.
- The HERMLE **SIEVA 3** is non explosion-proof and must therefore not be operated in explosion-endangered areas or locations. Centrifugation of flammable, explosive, radioactive, or such substances, which chemically react with high energy, is strictly prohibited. The final decision on the risks associated with the use of such substances is the responsibility of the user of the centrifuge.
- Never spin toxic or pathogenic material without adequate safety precautions, i.e. centrifugation of buckets / tubes without or with defective hermetic sealings is strictly prohibited. The user is obliged to perform appropriate disinfection procedures in case dangerous substances have contaminated the centrifuge and or its accessories. When centrifuging infectious substances, always pay attention to the General Laboratory Precautions. If necessary, contact your safety officer!
- It is prohibited to run the centrifuge with rotors other than listed for this unit.
- Under no circumstances open the lid of the centrifuge while the rotor is still running or rotating with a speed of $> 2\text{m/s}$

The following rules must be strictly adhered to:

- Do not operate the centrifuge in case it is not installed correctly.
- Do not operate the centrifuge when dismantled (e.g. without housing).
- Do not run the centrifuge when mechanical or electrical assembly groups have been tampered with unauthorized persons.
- Do not use accessories such as rotors and buckets, which are not exclusively approved by HERMLE Labortechnik GmbH, except commercially available centrifuge tubes made of glass or plastic.
- Do not spin extremely corrosive substances, as they may cause material damages and impair mechanical resistance.
- Do not operate the centrifuge with rotors or buckets, which show any signs of corrosion or mechanical damage.

The manufacturer is responsible for safety and reliability of the centrifuge, only if:

- The unit is operated in accordance with this instruction manual.
- Modifications, repairs or other adjustments are performed by HERMLE-authorized personnel and the electrical installation of the related location corresponds to the IEC-regulations.

1.2 Description

The Sieva-3 is a filtration centrifuge with microprocessor control. All parameters are accessible via buttons and selected with the control field. All pre-selected and current values will be shown permanently on the LCD-display.

The set parameters remain stored after the centrifugation.

The lid is locked and unlocked by an electromagnetic lid lock.

The centrifuge is powered by a maintenance-free and silent induction motor.

1.3 Delivery Package

- 1 Centrifuge Sieva -3
- 1 Instruction Manual Sieva-3
- 1 Hex key
- 1 Allen key
- 1 Power cord


Rotor(s) / Accessories will be packaged either inside the centrifuge or separately.

1.4 Warranty

The centrifuge has been subjected to thorough testing and quality controls. In the unlikely case of any manufacturing faults occurring, the centrifuge and rotors are covered by warranty for a period of two years from date of delivery. This warranty becomes invalid in case of mishandling, damage and negligence and further in case of usage of inappropriate spare parts and / or accessories or unauthorized modification of the unit.

Technical modification rights are reserved by the manufacturer in respect to technical improvement!

1.5 Operating and Display Elements

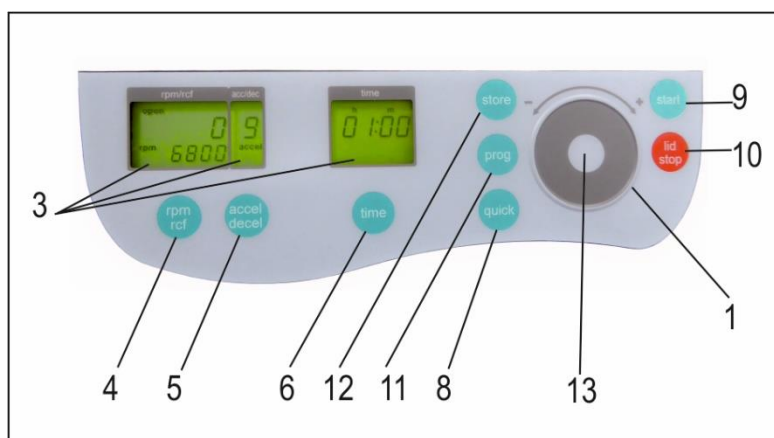

1	control field	Run Parameters
3	LCD	Control Panel Display
4	rpm/rcf	Speed / g-force
5	accel/decel	Acceleration / Deceleration
6	time	Centrifugation time
8	quick	Short Running
9	start	Start Centrifugation
10	lid/stop	Lid Release / Stop Centrifuge
11	prog	Retrieving Stored Programs
12	store	Program Store
13	LED light	Shows the status of the centrifuge


1.5.1 LED light

The LED light indicates the current operation state of the centrifuge. The following table shows all operating states.

Color of LED light	Operating state
Green	Run complete, lid is closed
Yellow	Lid opened, no sleep mode
Red flashing fast	Error message
Red flashing slow	Centrifuge runs
Yellow flashing slow	Sleep mode
Red – Green flashing slow	Standard settings menu

1.5.2 LCD-Display

The following picture shows the individual elements of the LCD-display.

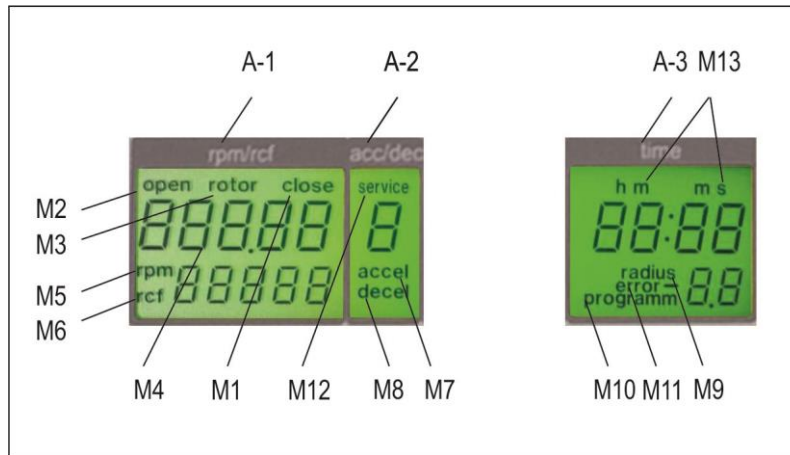

Display Fields:

- A-1 Display Field – "rpm/rcf"
- A-2 Display Field – "acc/dec"
- A-3 Display Field – "time"

Messages/Lettering of the Display Fields:

M1	"close"	M8	"decel"
M2	"open"	M9	"radius"
M3	"rotor"	M10	"program"
M4	Rotor-No.	M11	"error"
M5	"rpm"	M12	"service"
M6	"rcf"	M13	h m s
M7	"accel"		

1.6 Signs and Indications of the Centrifuge

1.6.1 General



Direction of Rotation – clockwise rotation for the rotor drive

1.6.2 Product Nameplate (Example)



Company Address: Hermle Labortechnik GmbH, Siemensstr. 25, D-78564 Wehingen

TYPE: Type Designation of the Product

REF: Order No. of the Product

SN: Serial Number of the Product



Date of Manufacture



Manufacturer



Instructions for Disposal, (see chapter 7, S. 26)

MAX. Drehzahl: Max. Allowed Speed of the Unit

KIN. EN.: Max. Kinetic Energy with Corresponding Rotor

U/I/f: Allowable Voltage / Max. Current / Frequency

P: Electrical Input Power



Operating Manual



RoHS-Compliance label (see chapter 7, S. 26)



Labeling identifies that all Standards and Guidelines are applied

1.6.3 Warning and Information Signs

Warning

Four carrier must be used at all times on four place swing out rotors or damage will occur to the centrifuge. Such damage will not be covered under the product warranty.

Four carrier must be used at all times on four place swing out rotors or damage will occur to the centrifuge. Such damage will not be covered under the product warranty.

Attention!!
Check the fastening
of the rotor nut before each run.
Achtung!!
Vor jedem Lauf Befestigungsschraube auf festen Sitz pruefen.

Attention! Check the fastening of the rotor nut before each run!



Vor manueller Notentriegelung
Netzstecker ziehen!
Disconnect main power plug
before manual emergency release!
Tirer fiche de contact avant une
déverrouillage d'arrêt d'urgence!

Take off mains plug before opening the housing or the emergency release



230 V/50-60 Hz

Power Input

2. INSTALLATION

2.1 Unpackaging

Model **Sieva-3** is supplied in a carton on a pallet. Remove the strap retainer and lift the carton. Now the centrifuge is standing free on the pallet. Lift the centrifuge on both sides with an appropriate number of helpers and place it on the laboratory table (see Figure 1).



Attention! Do not lift the centrifuge from under the lid or by the front panel!



Figure 1

The instruction manual must be kept with the centrifuge, at all times!

2.2 Space Requirements

The centrifuge should be installed on an even, solid surface, if possible on a laboratory cabinet / table or some other solid vibration free surface.

During centrifugation, the centrifuge must be placed in a way, that there is a minimum space of 30 cm/11.81in on each side of the unit, according to EN 61010-2-020 standards.

Do not place the centrifuge next to a window or a heater where it could be disposed to excessive heat, as the performance of the unit is based on an ambient temperature of 23°C/73.4°F.

2.3 Installation

Follow these steps:

- Check whether the power supply corresponds with the one specified on the manufacturer's rating label, mounted on the rear pane.
- The power connection for the centrifuge requires a **separate** one-site protection, with 16 A (Type K).
- In case of emergency, there must be an emergency switch off installed outside of the room, in order to disconnect the power supply from the unit.
- Connect the centrifuge, with the mains.
(The socket for the power cord must be easy to reach, respectively easy to disconnect).
- Switch on, by using the mains power switch (I).
- Open the lid, by using the button LID.
- Remove the transport securing device of the motor.

2.4 Installation of the rotor

2.4.1 Mounting of the centrifugal basket

Place the basket onto the shaft mounting. Make sure that the driving pins on the shaft mounting and the borings on the bottom of the basket are in the correct position. Hold the basket with one hand and secure the fixing nut with the help of the hex key by turning it clockwise (see Figure 2).



To ensure safe operation, the motor mounting must be regularly checked for tightness and secured, if necessary. Therefore remove the fixing nut from the basket. Hold the basket with one hand and tighten the motor mounting with the help of allen key by turning it clockwise (see Figure 3).



Figure 2



Figure 3

2.4.2 Überladung von Rotoren

The maximum load permitted for the rotor is determined by the manufacturer, as well as the maximum speed allowed for this rotor (see label on rotor), must not be exceeded.

The liquid the rotors are loaded with, should have a max. homogeneous density of 1.2 g/ml or less when the rotor is running at maximum speed

In order to spin liquids with a higher density, the speed has to be reduced, according to the following formula:

$$\text{Reduced speed } n_{\text{red}} = \sqrt{\frac{1,2}{\text{higher density}}} \times \text{max. speed } (n_{\text{max}}) \text{ of the rotor}$$

Example:

$$n_{\text{red}} = \sqrt{\frac{1,2}{1,7}} \times 4.000 = 3.360 \text{ rpm}$$

2.4.3 Removing of the rotor

Hold the basket with one hand and untighten the fixing nut of the basket by turning the hex key counter clockwise. Lift the rotor from the motor mounting.

2.4.4 Functional principle of the baskets

The baskets of Sieva-3 are available in two versions. As a perforated and a non-perforated basket.

The perforated basket 221.07 V02 is used for the filtration method. Therefore, a filter with desired pore size is placed into the inner wall of the basket. Due to the centrifugal force, small particles of the solution remain suspended in the filter and the filtered liquid is pressed through the holes of the basket (see Figure 4).

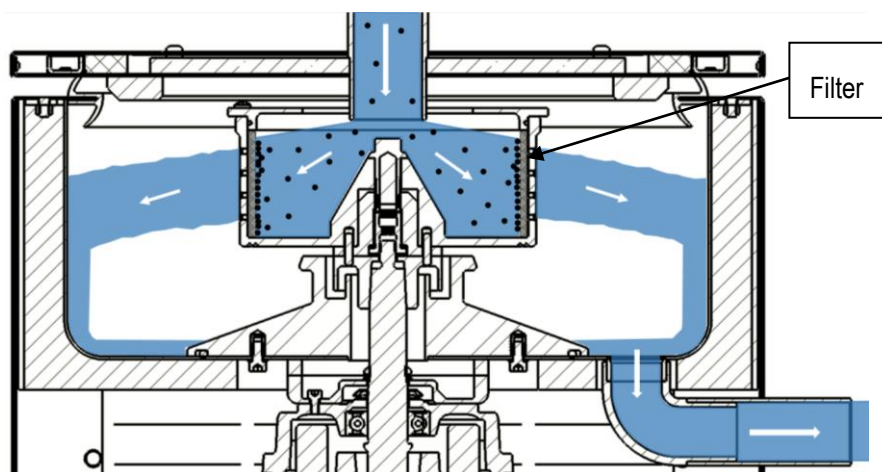


Figure 4: Filtration method

The non-perforated basket 221.07 V04 is used for decantation method. After adding of the solution, the sediment is pressed against the inner wall of the basket and stands upright due to the centrifugal force. The light liquid is discharged via the opening in the lid of the basket (see Figure 5).

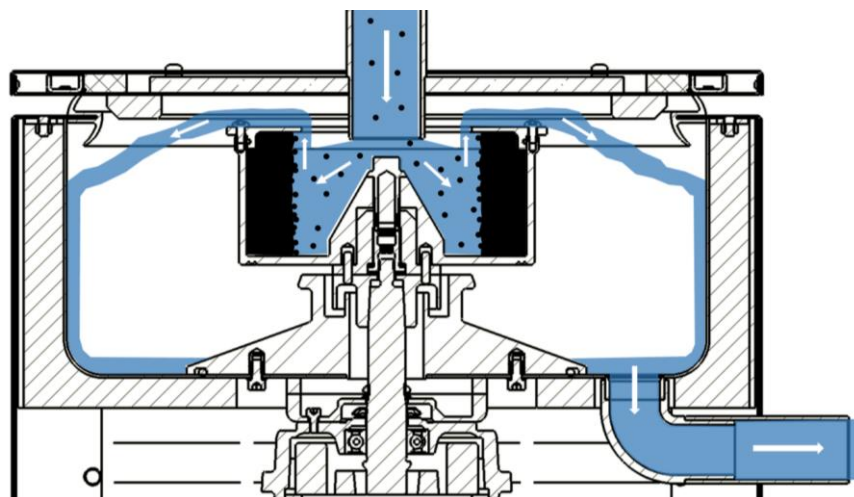


Figure 5: Decantation method

3. BASIC ADJUSTMENTS

3.1 Adjustment of the Rotor Type

In the centrifuge model **SIEVA-3** no rotor adjustment can be done.

3.2 Access to the Mode: “Standard settings“

While starting this model, the following basic settings can be performed:

- Acoustic Signal Turn On / Off
- Volume Pre-Selection of Sound Signal
- Song Selection of Sound Signal "end of run"
- Keyboard Sound On / Off
- Sleep Mode On / Off

The following operating data can be retrieved in this mode:

- Number of Starts
- Operating Hours of Centrifuge
- Operating Hours of Motor
- Software Version
- Frequency Converter Software
- Error List
- Function of the Imbalance Switch
- Operation of Keyboard
- Hardware Version
- Serial Number and year of manufacturing of the control board
- Update touch control panel

Open the centrifuge lid and shut off the main switch. Now switch on again the main switch. For approximately 3 seconds “SIEVA“ will be shown in the display. Press during this time the keys “time“ (6) and “lid“ (10) simultaneously. As a result, a display test is administered for approx. 3 seconds. All possible indications will appear at the same time (see Figure 6). The LED light is flashing alternately with green and red color.

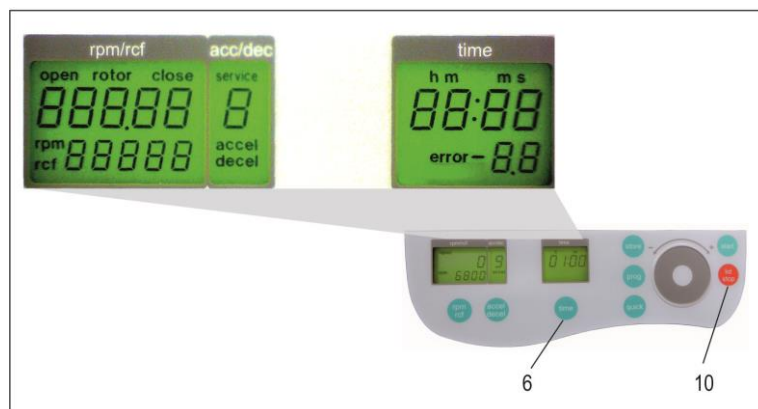


Figure 6



ATTENTION:

- The normal program mode can be changed back again by switching off the centrifuge, for a short period!
- All changed settings must be confirmed by the key, “start” (9). A confirmation screen will appear with the word, “store”, in the display “rpm/rcf” (A-1) - Only then the pre-selections are valid!

3.3 Acoustic Signal Turn On / Off

Proceed as illustrated, under point 3.2, to enter this program mode, press the key; "accel/decel" (5). In the display, "accel/decel" (A-2) flashes the word, "service". Select the letter, "L" with the control field (1). As a result, appearing in the display "rpm/rcf" (4), are the words, "On Sound". By pressing the key, "rpm/rcf" (4), the word "On" flashes, and the sound can be switched off with the control field (1), (see Figure 7).

After the settings have been stored by user, the normal program mode can be changed back again by switching off the centrifuge, for a short period.

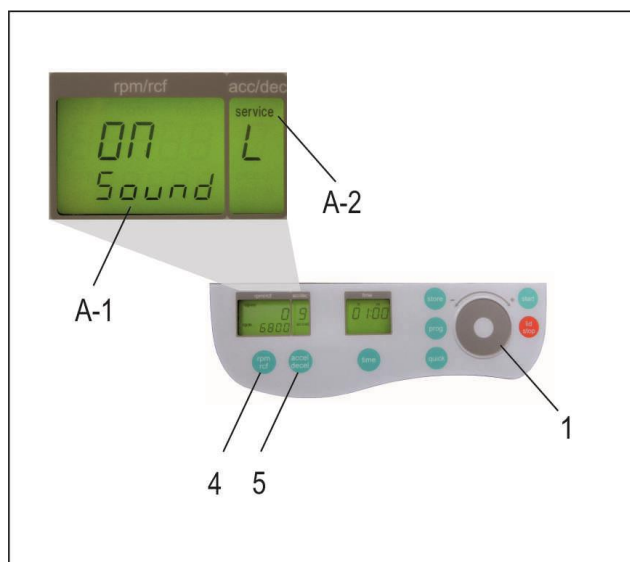


Figure 7

3.4 Volume Pre-Selection of Sound Signal

Proceed as illustrated, under point 3.2, to enter this program mode, press the key "accel/decel" (5). In the display, "accel/decel" (A-2), flashes the word, "service". Select the letter, "U" with the control field (1). As a result, appearing in the display, "rpm/rcf" (A-1), are the words, "Vol=0- 9/Sound". By pressing the key, "rpm/rcf" (4), the desired volume can be adjusted between 0 (low) and 9 (loud), with the adjustable knob (1), (see Figure 8).

After the settings have been stored by user, the normal program mode can be changed back again by switching off the centrifuge, for a short period.

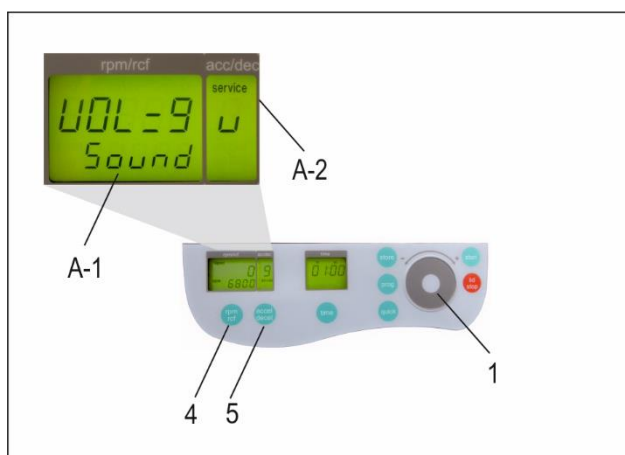


Figure 8

3.5 Song Selection – End of Run

Proceed as illustrated, under point 3.2, to enter this program mode, press the key, "accel/decel" (5). In the display, "accel/decel" (A-2) flashes the word, "service". Select the letter, "G" with the control field (1). As a result, appearing in the display, "rpm/rcf" (A-1), the word " SonGo/Sound". After pressing the key „rpm/rcf“ (4), select a song with the adjusting knob (1), (see Figure 9).

After the settings have been stored by user, the normal program mode can be changed back again by switching off the centrifuge, for a short period.

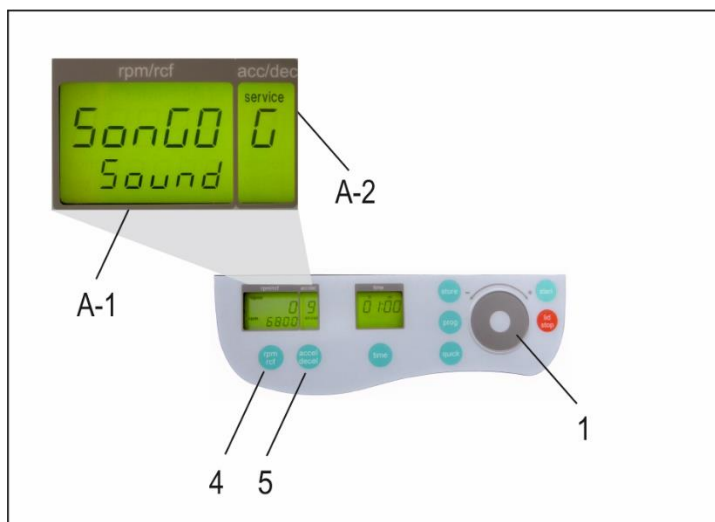


Figure 9

3.6 Keyboard Sound Turn On / Off

Proceed as illustrated, under point 3.2, to enter this program mode, press the key, "accel/decel" (5). In the display; "accel/decel" (A-2) flashes the word, "service". Select the letter, "b" with the control field (1). As a result, appearing in the display, "rpm/rcf" (A-1), the word "ON/BEEP". By pressing the key, "rpm/rcf" (4), the keyboard sound (On) or (Off) can be turned on, with the adjustable knob (1), (see Figure 10).

After the settings have been stored by user, the normal program mode can be changed back again by switching off the centrifuge, for a short period.

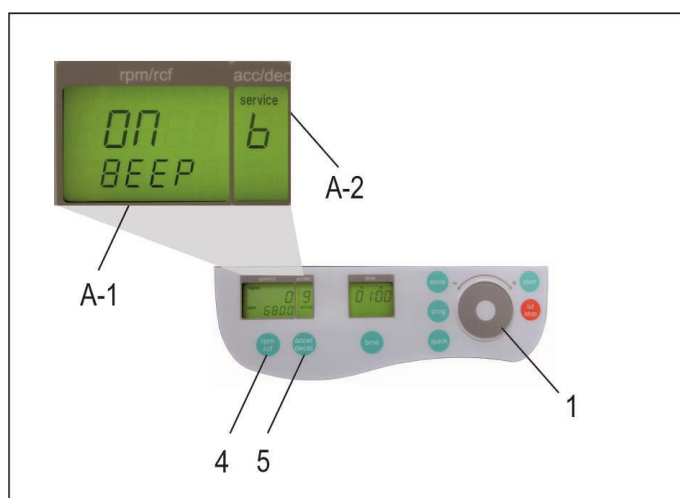


Figure 10

3.7 Sleep Mode Setting

After not using the centrifuge for 5 minutes, the centrifuge automatically falls into sleep mode. The display is switching off and the LED light is flashing slowly with yellow color. By pressing any key, the centrifuge will be reactivated. The sleep mode can be deactivated or set in the range between 1 – 60 minutes.

Proceed as illustrated, under point 3.2, to enter this program mode, press the key, "accel/decel" (5). In the display, "accel/decel" (A-2) flashes the word, "service". Select the letter "I" with the control field (1). As a result, appearing in the display, "rpm/rcf" (A-1), the word "SLEEP". After pressing the key „rpm/rcf“ the sleep mode can be turned off or pre-set to the needed value by turning the control field (1).

After the settings have been stored by user, the normal program mode can be changed back again by switching off the centrifuge, for a short period.

3.8 Retrieving Operation Data

In the function, "Basic Adjustments" the operating data, of the centrifuge, can be retrieved. Please proceed as described, under point 3.2, to enter this program mode.

Press the key "accel/decel" (5). In the display, "accel/decel" (A-2) flashes the word, "service".

With the control field (1) the following information can be retrieved:

A	= Previous Starts of the Centrifuge
H	= Previous Operating Hours
h	= Running Time of the Motor
S	= Software Version
r	= Frequency Converter Software
E	= List of Previous Error Messages
F	= Function of the Imbalance Sensor
P	= Operating of Keyboard
d	= Hardware-Version
y	= Serial Number and year of manufacturing of the control board
n	= Update touch control panel (only by trainer service staff)

The list of the last 99 error messages can be looked over by pressing the key "rpm/rcf" (4) and leaf through it with the control field (1). The respective error codes appear in the display "rpm/rc" (A-1) (see Figure 11). The first two numbers indicate the appeared errors ongoing from 00 to 99, the last two numbers indicate the error code. Please refer to, "Table 4: " (see APPENDIX p. V)

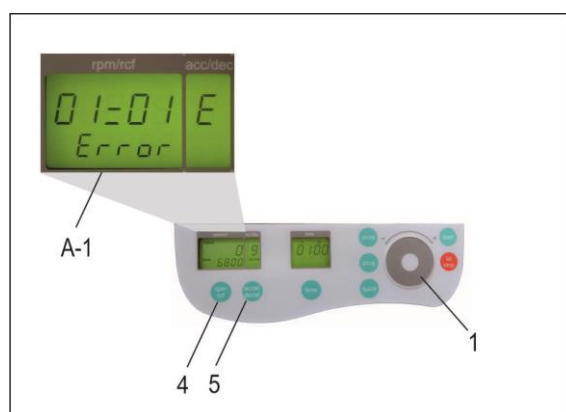


Figure 11

4. OPERATION

4.1 Power Switch

The power switch is located on the left site of the centrifuge housing (see Figure 12).



Figure 12: Power Switch



ATTENTION:

After turning on the power switch, open the centrifuge lid first, before starting the centrifuge.

4.2 Centrifuge Lid

4.2.1 Lid Release

After the run, properly close the lid of the centrifuge, appearing in the display, "rpm/rcf"(A-1) with the word, "close" (M1). At the same time the actual rotor type, for example "221.57", is shown in the display. By pressing the key, "lid" (10), the lid of centrifuge can be released. As soon as the lid is completely released, the word, "open" (M2) appears. The lid of the centrifuge is now able to be opened.

For all number marked text, please refer to Figure 13.

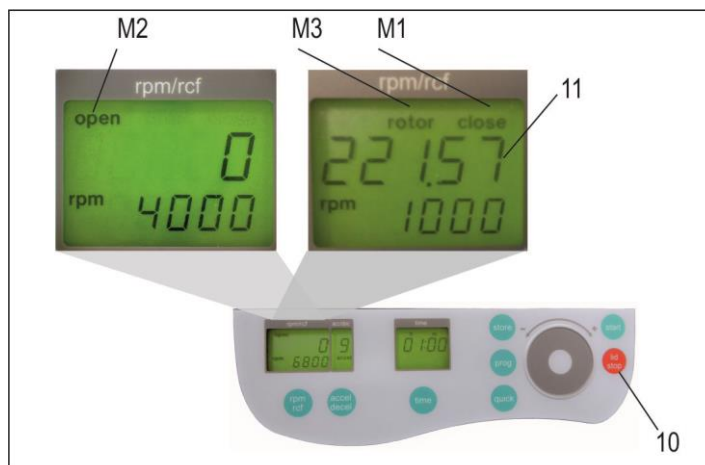


Figure 13

4.2.2 Lid Lock

The lid must only be closed slightly. After closing the lid, the word "open" (M2) will no longer be displayed. As a sign that the centrifuge is ready for starting, appearing in the display, "rpm/rcf" (A-1), the word "close" (M1). Simultaneously, the word "rotor" (M3) appears, as well as the code number of the rotor, which is in the centrifuge, along with all rotor specific data, for example: max. speed, acceleration etc., are available.

For all number marked text, please refer to Figure 13.



ATTENTION:

Before closing the lid, please check if the rotor is tightened.

4.3 Pre-Selection

4.3.1 Pre-Selection of Speed and RCF-Value

By selecting the key, "rpm/rcf" (4), pre-selection is activated. By pressing the key once, the word "rpm" (M5) flashes. By pressing the key again, the pre-selection of the centrifugal forces can be chosen. The flashing word, "rcf" (M6), will appear.

The desired values can be selected, with the control field (1). In the display (A-1), the regulated value is shown permanently: before, during and after the run.

This pre-selected value will be stored, as long as a new pre-selection is made.

For all number marked text, please refer to Figure 14.

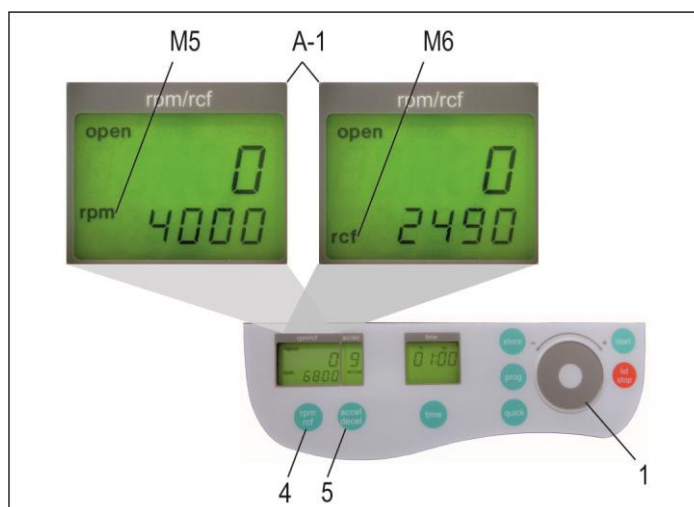


Figure 14

As long as no rotor is inserted, the speed is adjustable between 200 rpm and the maximum revolution of the centrifuge.

If there is a rotor in the centrifuge, the speed can only be pre-selected up to the maximum permissible revolution of that rotor. It is the same with the pre-selection of the RCF-Value. The setting range is between 20 xg and the maximum permissible centrifugal force of the rotor.

The permissible max. speed of the SIEVA-3 with the rotor 221.07 V02/04 is **10000 rpm** or **7825 xg**.

4.3.2 Pre-Selection of Running Time

The running time can be pre-selected in 3 different ranges: from 10 seconds up to 99 hours 59 minutes.

1. Range from: 10 seconds up to 59 minutes 50 seconds, in steps of 10 seconds
 2. Range from: 1 hour up to 99 hours 59 minutes, in steps of 1 minutes
 3. Range: Continuous Run "cont", can be interrupted by the key, "stop" (10).
- The running time can be pre-selected, with the lid opened or closed.
 - To activate the setting of the running time, press the key "time" (6).
 - In the display, "time" (A-3) flashes the indication: "m : s" or "h : m", depending on the previous setting.

To set the desired value, use the control field (1). After exceeding 59 min 50 sec, the indication changes automatically to, "h : m". After exceeding 99 hours 59 min, the word "cont" appears in the display, "time" (A-3).

The continuous run can only be interrupted by pressing the key, "stop" (10). The time counts down, as soon as the set speed is reached.

For all number marked text, please refer to Figure 15.

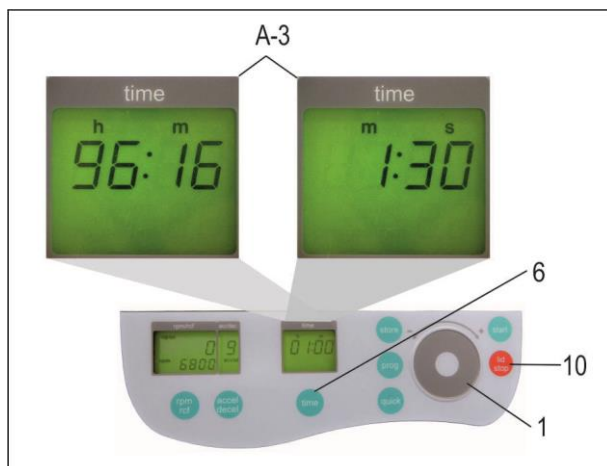


Figure 15

4.3.3 Pre-Selection of Brake Intensity and Acceleration

Selecting the key, "accel/decel" (5), this function is activated.

By pressing the key once, the word "accel" (M7) flashes, in the display "accel/decel" (A-2). The desired acceleration can be pre-selected, with the control field (1). The value 0 is equivalent to the lowest acceleration and the value 9 is equivalent to the highest acceleration.

By pressing the key "accel/decel" (5) twice, in the display "accel/decel" (A-2), indicates the word "decel" (M8). Now the desired brake intensity can be pre-selected, with the control field (1). The value 9 is equivalent to the shortest possible brake time and the value 0 to longest possible brake time.

For all number marked text, please refer to Figure 16.

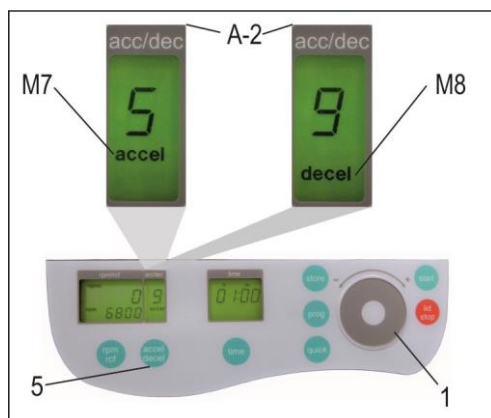


Figure 16

This table shows the acceleration and deceleration stages 0 and 9 of the empty baskets:

Rotor-Number	Acceleration		Deceleration	
	Stage 0	Stage 9	Stage 0	Stage 9
221.07 V02	135	15	450	18
221.07 V04	135	15	640	18
in seconds				
Acceleration time from 0 min ⁻¹ - n _{max}			Deceleration time from n _{max} - 0 min ⁻¹	

4.4 Program

4.4.1 Storage of Programs

The program stores up to 99 runs, with all relevant parameters, including the used rotors. Any free program number is available and can be retrieved.

Put the desired rotor into the centrifuge. By pressing the key, "prog" (11), in the display "time" (A-3) appears the word "program", (see Figure 17). With the adjustable knob (1), choose the desired program number.

If a program number is already occupied, in the display "rpm/rcf" (A-1), the words "rotor" (M3) and e.g. "221.57" (M4) will appear, (see Figure 17). Free program numbers will appear as 0.

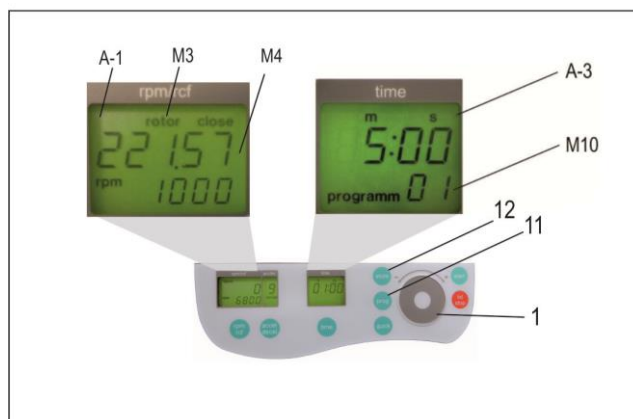


Figure 17

Close the lid of the centrifuge, now proceed as described above, to set all important run parameters. If the lid isn't closed, when storing the program in the display "rpm/rcf" (A-1), flashes alternately the word "FirSt" and "CLOSE Lid" (see Figure 18). When starting the run without storing the program, in the display "rpm/rcf" (A-1), flashes alternately the word "First" and "PrESS StoreE", (see Figure 19).



Figure 18



Figure 19

For alteration of data, press the key "store" (12), for approx. 1 second. If the program is stored correctly, the word "StorE" appears in the display "rpm/rcf" (A-1). As a result, the word "program" (M10) disappears. As soon as the key "store" (12) is no longer displayed, the word "programm xx" (M10) reappears, (the xx stands for the chosen program place).

If all program numbers are occupied, take an old number that is not needed any longer and replace it with the new parameters.

4.4.2 Recall of Stored Programs

To recall stored programs, press the key "prog" (11), with the lid already closed. Inside the display "time" (A-3), appears "programm --"(M10). With the control field (1), pre-select the desired program number. In the respective displays, the stored values, for that program, will appear.

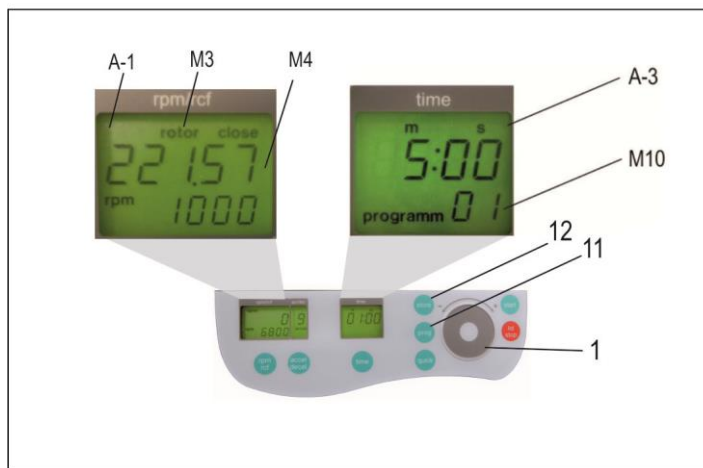


Figure 20

4.4.3 Leaving Program Mode

To leave the program mode, press the key, "prog" (11). Then, inside the display "time", appears the word "programm".

Set the display to "programm--" (M10) with the control field (1).

For all number marked text, please refer to Figure 20.

4.5 Starting and Stopping the Centrifuge

4.5.1 Starting the Centrifuge

Start the centrifuge with either the "start" key (9), or the "quick" key (8). With the "start" key (9), stored runs or runs with manually pre-selected parameters can be started. During the centrifugation the LED light is flashing slowly with red color. When the respective pre-selected running time has ended, the centrifuge will stop automatically.

By pressing the "quick" key (8), the centrifuge accelerates up to the pre-selected revolution. In the display "time" (A-3), the passed running time is indicated from the moment the "quick" key (8) is pressed. By releasing the "quick" key (8), the centrifuge stops and the running time is indicated, until the lid is opened.

For all number marked text, please refer to Figure 21.

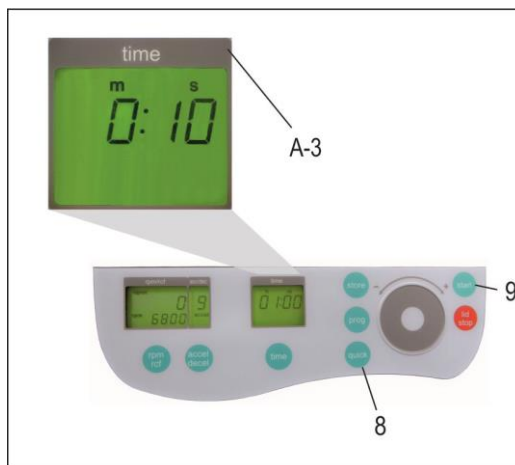


Figure 21

4.5.2 Stopping the Centrifuge

With the "stop" key (10), the run time can be interrupted, at any time, (see Figure 22). After pressing the key, the centrifuge decelerates with the respective pre-selected intensity, down to a standstill. After the centrifugation is finished, the LED lights with green color.



Figure 22

4.6 Imbalance Detection

With fast filling of the basket in the acceleration phase, an imbalance can occur. In this case, the centrifuge switches off the drive immediately and the basket decelerates to standstill.

The display "time" (A-3) will show the word "error" (M11) and the number "01" appears.

When inside the display "time" (A-3), the word "error" along with the number "02" appear, (see Figure 23) the cause may be as follows. -> The imbalance switch is defective.

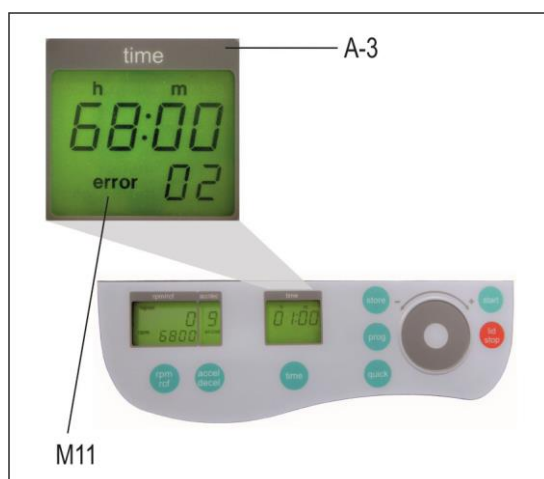


Figure 23

4.7 Volume Flow



The volume flow of the medium to be analysed may be max. 3.5 l/min. When the flow rate is higher, the temperature of the drive increases until it is switched off automatically for safety reasons.

4.8 External Temperature Control

The Sieva-3, with the preparation for external temperature equalization, is equipped with a copper pipe winding around the centrifuge chamber. On the side of the housing there are screw connections with an internal thread of 1/4". Hence, nipples of various sized can be used.



The tempering liquid may have a max. temperature of 90°C. The liquid may be pumped through the copper pipe winding with a max. pressure of 5 bar.

5. MAINTENANCE

5.1 Maintenance and Cleaning

General Care:

Maintenance of the centrifuge is dependent on prolonging the life of the rotor, the rotor chamber and the rotor accessories. Please be sure to clean the accessories, especially the sealing of the aerosol-tight rotors and insert bolts, of swing out rotors. Following, lubricate the bolts or sealing, with the recommended HERMLE Rotorgrease - Order No.: 38-5656.

Please pay special attention to anodized aluminum parts. Breakage of rotors can be caused by the slightest damages.

In case of rotors, buckets or tube racks becoming in touch with corrosive substances, the affected area must be cleaned, thoroughly.

Corrosive substances, such as, must be avoided: alkalis, alkaline soap solutions, alkaline amines, concentrated acids, solutions containing heavy metals, water-free chlorinated solvents, saline solutions, e.g. salt water, phenol, halogenated hydrocarbons.



Cleaning Units, Rotors, Accessories:

- Turn the device off and disconnect from the power supply, before beginning any cleaning or disinfecting. Do not pour liquids into the housing interior.
- Spray disinfectant on the device.
- Thorough cleaning not only has its purpose in hygiene, but also in avoiding pollution based corrosion.
- In order to avoid damaging anodized parts, such as rotors, reduction plates etc.; only pH-neutral Detergents, with a pH-value of 6-8, may be used for cleaning. Alkaline cleaning agents must not be used, (pH-value > 8).
- After cleaning, please ensure all parts are dried thoroughly, either by hand or in a hot-air cabinet (Max. Temperature + 50°C/122°F).
- It is necessary to coat anodized aluminium parts with anti-corrosion oil regularly, in order to increase their life-span and reduce corrosion predisposition.
- Due to humidity or not hermetically sealed samples, condensation may form. The condensation has to be removed from the rotor chamber, with a soft cloth regularly



The maintenance procedure has to be repeated every 10 to 15 runs, but at least once a week!

- Connect the unit to the power supply only after the equipment is completely dry.
- Do not implement disinfection with UV-, beta- and gamma-rays or other high energy radiation.
- Metal rotors can be autoclaved.
- Rotor lid can also be autoclaved, (Max. 121°C/250°F, 20 min).
- The tube racks are made of PP and **cannot** be autoclaved, at 134°C/273°F.

Cleaning and Disinfecting of the Unit

1. Open the lid, before turning off the unit. Disconnect from the power supply.
2. Open the rotor nut, by turning the rotor key counter-clockwise.
3. Remove the rotor from the motor mounting
4. For cleaning and disinfection of the unit and the rotor chamber, use the above mentioned cleaner.
5. Clean all accessible areas of the device and its accessories, including the power cord, with a damp cloth.
6. Wash the rubber seals and rotor chamber thoroughly, with water.
7. Rub the dry rubber seals with glycerol or talc, to prevent these from becoming brittle. Other components of the unit, e.g. the motor shaft or rotor, should **not** be greased.
8. Dry the motor shaft with a soft, dry and lint-free cloth.
9. Examine the unit and accessories for damage.

Remove adherent dust, at least every 6 months, from the ventilation slots in the centrifuge, by using a soft brush.

Cleaning and Disinfection of the Rotor

1. Clean and disinfect: the rotors, rotor lids and adapters, with the cleaner previously mentioned above.
2. Use a bottle brush to clean and disinfect the rotor bores.
3. Rinse the rotors, rotor lid and adapter, with clear water. Particularly, the drillings of the angle rotors.
4. When drying the rotors and accessories, set on a towel. Place the angle rotors, with bores down, to dry.
5. Dry the rotor cone with a soft, dry and lint-free cloth, check for damage. Do not grease the rotor cone.


Disinfection of Aluminum Rotors

In case of infectious material spilling into the centrifuge, the rotor and rotor chamber have to be disinfected, promptly after the run. Rotors may be autoclaved, at a maximum temperature of 121°C/250°F.

Disinfection of PP-Rotors

Autoclaving

The recommended time for autoclaving: 15 – 20 min at 121°C/250°F, (1 bar)


 **ATTENTION: The sterilization time of 20 min. must not be exceeded. Continuous sterilization will cause reduction in the mechanical resistance, of the plastic material.**

Before autoclaving the PP-rotor and adapter, thoroughly clean to avoid the burning of dirty residue.

Please disregard any consequences of chemical residues to plastic materials, at ambient temperatures. At high temperatures, autoclaving residue may corrode and destroy the plastic. The objects must be thoroughly washed with distilled water, after the cleaning, but before the autoclaving. Residues of any cleaning liquids, may cause fissures, whitening and stains.


Gas Sterilization

Adapters, bottles and rotors may be gas sterilized, with Ethylenoxyd. According to the duration of the application, allow items to properly air out, after the sterilization and before usage

 **ATTENTION: The temperature may rise during the sterilization; rotors, adapters and bottles should not be fully closed, keep completely unscrewed.**

Chemical Sterilization


Bottles, adapters and rotors may be treated, with the usual liquid disinfectants.

 **ATTENTION: Before applying any other, Cleaning Resp. Decontamination Method, other than what was recommended by the manufacturer, contact the manufacturer to ensure that it will not damage the unit or the rotor.**

Glass Breakage

With high g-values, the rate of glass tube breakage increases. Glass splinters have to be removed immediately from rotor, buckets, adapters and the rotor chamber itself. Fine glass splinters will scratch and therefore damage the protective surface coating of a rotor. If glass splinters remain in the rotor chamber, fine metal dust will build up, due to air circulation. This very fine, black metal dust will severely pollute the rotor chamber, the rotor, the buckets, and the samples.

If necessary, replace the adapters, tubes and accessories, to avoid further damage. Check the rotor bores regularly, for residue and damage.

 **ATTENTION: Please check the relevant specifications of the tubes centrifuges with the manufacturer!**

Lifetime of Rotors, Round and Rectangular Buckets, Accessories

Rotors and rotor lid made of aluminum or stainless steel, have a operating time of **max. 7 years** from first use.

Transparent rotor lids and caps, made of PC or PP, as well as rotors, tube racks and adapters of PP, have a maximum operating time of up to **3 years**, from first time use.

Conditions For the Operating Time:

Proper use, damage-free condition, recommended care.

6. TROUBLE SHOOTING

6.1 Error Messages: Problem / Solution

The error messages are listed to help localize possible errors faster.

The possible error referred to in this chapter may not always be the case, as they are only theoretically occurring errors and solutions.

Always keep us informed about any kind of error occurring, which is not listed in this chapter. With this information provided, we are able to improve and complete this operation manual.

Many thanks in advance for your support.

HERMLE Labortechnik GmbH

6.2 Survey of Possible Error Messages and Solutions

6.2.1 Lid Release during Power Failure (Emergency Lid Release)

In case of power failure or malfunction, the lid of the centrifuge can be opened manually, in order to protect samples.

Please proceed as follows:



- **Switch the centrifuge off and unplug the power cord, wait until the rotor is at a standstill, (this may take several minutes).**
At the right side of the centrifuge, there is a plastic stopper, (Figure 24 Remove this stopper, which is connected to the lid lock, horizontally from the housing until the centrifuge lid opens.
- Now, open the lid of the centrifuge.



Figure 24

6.2.2 Description of the Error Message System

The error message, "error" (M11), is shown in the "time" (A-3) display, (see Figure 25). For more detailed information, refer to Table 4: , (see Appendix P.V).

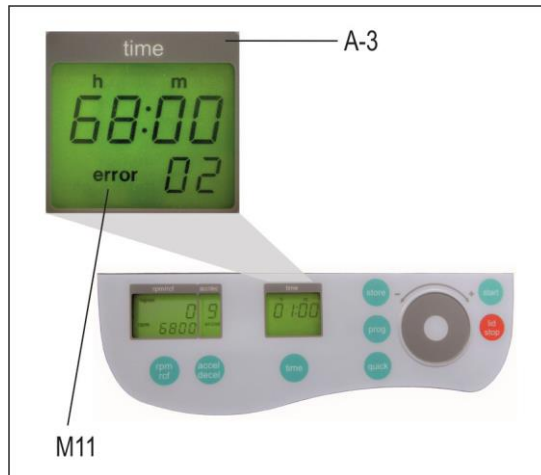


Figure 25

6.2.3 Procedure while error 14

If Error 14 occurs, there is a problem with the speed sensor. The centrifuge lid is closed for undefined period of time and in the "rpm/rcf" (A-1) display shows the lettering "USEr GuidE" (see Figure 26).

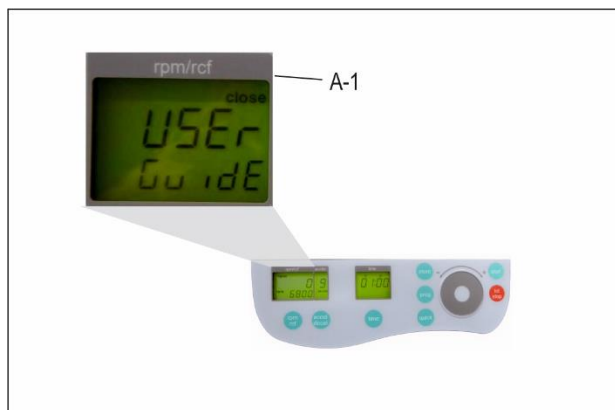


Figure 26



To reopen the centrifuge lid, switch off the device and wait until the rotor has come to a **standstill!** Take from chapter 4.3.3 the maximum deceleration time of the respective rotor. Level 0 corresponds to unbroken rundown, which occurs at error 14. If the centrifuge lid is opened before standstill of the rotor, a following error can occur.

Once the rotor has come to a standstill, open the centrifuge lid with the emergency release. Proceed as described in chapter 6.2.1, P. 24. After opening the centrifuge lid, switch on the device again. Error 14 and the lettering "USEr GuidE" should be eliminated.

If this error occurs again, please contact the service.

7. TRANSPORT, STORAGE AND DISPOSAL

Transport

- Before transporting, take out the rotor.
- Only transport the unit in its' original packaging.
- Use a transport aid, for transporting over longer distances, to fix the motor shaft.

	Air Temperature	Rel. Humidity	Air Pressure
General Transportation	-25 to 60 °C	10 to 75 %	30 to 106 kPa
	-13°F to 140°F		

Storage

During storage of the centrifuge, the following environmental conditions must be observed:

	Air Temperature	Rel. Humidity	Air Pressure
Transport Packaging	-25 to 55 °C	10 to 75 %	70 to 106 kPa
	-13°F to 131°F		

Disposal

In the event of disposing of the product, please observe the applicable legal requirements.

Information on the disposal of the electrical and electronic devices, in the European Community:

disposal of the electrical devices is regulated within the European Community, by national regulations, based on EU Directive 2002/96/EC pertaining to waste electrical and electronic equipment (WEEE).

In accordance with this, any devices delivered after 08/13/2005 on a business-to-business basis, including the product, may no longer be disposed of, as a household waste. To document this, the devices have been marked with the following identification:



Because disposal regulations may differ from one country to another, within the EU, please contact your supplier, if necessary.

RoHS II Compliance

HERMLE Labortechnik GmbH, Siemensstraße 25, 78564 Wehingen, hereby declares and certifies that all components manufactured are RoHS II compliant, according to the definition and restrictions given by the European Parliament Directive 2011/65/EC. This restricts the use of certain hazardous substances in electrical and electronic equipment.

8. APPENDIX

EC – Conformity Declaration	II
Tabelle 1: Technical Data.....	III
Table 2: Max. Speed and RCF-Values for Permissible Rotors	IV
Table 3: Permissible Net Weight	IV
Table 4: Error Messages	V
Table 5: Abbreviations	V

EC – Conformity Declaration

**EG Konformitätserklärung
EC Conformity Declaration****Hermle Labortechnik GmbH - Siemensstr. 25 - D-78564 Wehingen – Germany**

Das bezeichnete Produkt entspricht den einschlägigen grundlegenden Anforderungen der aufgeführten EG-Richtlinien und Normen.
Bei einer nicht mit uns abgestimmten Änderung des Produktes oder einer nicht bestimmungsgemäßen Anwendung verliert diese
Erklärung ihre Gültigkeit.

The Product named below fulfills the relevant fundamental requirements of the EC directives and standards
listed. In the case of unauthorized modifications to the product or an unintended use this declaration
becomes invalid.

Produkttyp
Product type

Laborzentrifugen mit Zubehör (sonstige Produkte)
Laboratory centrifuge with accessories (other device)

Typenbezeichnung
Typ designation

SIEVA-3

Einschlägige EG-Richtlinien / Normen
Relevant EC directives / standards

**2006/95/EG, 2004/108/EG, RoHS II 2011/65/EG,
EN 61 010-1; EN 61 010-2-020; ISO 9001:2008**

Wehingen, den 09.09.2018

(gültig bis/valid until 21.12.2018)

HERMLE
LABORTECHNIK 



Alexander Hermle
Geschäftsführer - Managing Director

Tabelle 1: Technical Data

Manufacturer		HERMLE Labortechnik GmbH 78564 Wehingen			
Model Type		Sieva-3			
Dimension					
Width		38 cm			
Depth		50 cm			
Height		32 cm			
Weight without rotor		37 kg			
max. Speed		10000 rpm			
max. Volume		500 ml			
max. RCF		7825 x g			
Allowed density		1,2 kg/dm ³			
Allowed kinetic energy		6300 Nm			
Mains Power Connection AC		230 V / 50 Hz 1 ph		120 V / 60 Hz 1 ph	
Voltage Fluctuation		± 10%		± 10%	
Current Consumption		3,5 A		6,8 A	
Power Consumption		0,8 KW		0,81 KW	
Radio Interference		IEC 61326-1			
Audit Requirement (BGR 500)		no			
Ambient Conditions (EN/IEC 61010-1)					
- Environment - High - Ambient Temperature - Max. Relative Humidity - Overvoltage Category (IEC 60364-4-443) - Degree of Contamination		for indoor use only Use up to an altitude of 2000 m above MSL 2°C up to 35°C Max. relative humidity 80 % for temperatures up to 31°C, decreasing linearly to 50 % relative humidity up to 35°C. II 2			
Class of Protection I		Class of protection (DIN EN 60529) IP 20			
Not suitable for use in hazardous environments					
EMC		EN/IEC	FCC Class B	EN/IEC	FCC Class B
Interference Emission		61326-1		61326-1	
		Class B		Class B	
Noise Level (depending on the rotor)		64-70 dB(A), depending on the rotor and the filter paper			
Write from the Operator					
Inventory-No.:					
Monitoring-No.:					
Environment:					
Maintenance Contract:					
Responsible Service Office			HERMLE Labortechnik GmbH		Dealer or Service Office
			Siemensstraße 25		
			78564 Wehingen		
			Tel.: (49)7426 / 96 22-17		
			Fax: (49)7426 / 96 22-49		
Responsible Dealer					

Table 2: Max. Speed and RCF-Values for Permissible Rotors

Rotor Number	Max. Speed	RCF Value
221.07 V02	10000 rpm	7825 xg
221.07 V04	10000 rpm	7825 xg

Table 3: Permissible Net Weight

Rotor Number	Max. Speed	Permissible Net Weight
221.07 V02	10000 rpm	600 g
221.07 V04	10000 rpm	600 g

Table 4: Error Messages

Error-No.:	Description
1	Imbalance arose
2	Imbalance sensor is defective
4	Imbalance switch has been activated for longer than 5 seconds
8	Transponder in the rotor is defective
14 USEr GuidE	Leap of speed is too large between two measurements. Centrifuge lid is closed for undefined period of time
16	Standstill detection defective
33	Open lid while motor is running
40	Communication with frequency converter disturbed during start
41	Communication with frequency converter disturbed during stop
42	Short circuit in the frequency converter
43	Under-voltage frequency converter
44	Overvoltage frequency converter
45	Over temperature frequency converter
46	Over temperature motor
47	Over current frequency converter
48	Timeout between control unit and frequency converter
49	Other error frequency converter
55	Over Speed
99	Rotor not allowed in this centrifuge
FALSE	Inserted rotor does not exist in the program
rotor no	Rotor is not detected

Table 5: Abbreviations

Symbol / Abbreviation	Unit	Description
U (=rpm)	[min ⁻¹]	Revolutions per Minute
RZB(=rcf)	[x g]	Relative Centrifugal Force
PP	-	Polypropylene
PC	-	Polycarbonate
accel	-	Acceleration
decel	-	Deceleration
prog	-	Program



Redemption Form / Decontamination Certificate

Decontamination Certificate of Goods Returned upon Delivery

Enclose all returned shipping items and modules necessary!

The complete, full declaration about the decontamination is a prerequisite for the assumption and further processing of the return. If no corresponding explanation is enclosed, we carry out decontamination with costs at your expense.

Surname; Last Name: _____
Organization / Company: _____
Street: _____
ZIP CODE: _____ **place:** _____
Telephone: _____ **fax:** _____
E-Mail: _____

Please fill out in block
capitals!

Pos.	Quantity	Decontaminated Object	Serial No.	Description/Comment
1				
2				
3				
4				

Are the parts listed above in touch with the following substances?

Health endangering watery solutions, buffers, acids, alkalis:..... ☐ Yes ☐ No

Potentially infectious agents: ☐ Yes ☐ No

Organic reagents and solvent: ☐ Yes ☐ No

Radioactive substances: ☐ α .. ☐ β .. ☐ γ .. ☐ Yes ☐ No

Health endangering proteins: ☐ Yes ☐ No

DNA: ☐ Yes ☐ No

Have these substances reached the equipment/assembly? ☐ Yes ☐ No

If so, which ones:

Description of the measures for the decontamination of the listed parts:

I confirm the proper decontamination:

Company/Dept . _____ Place and Date: _____

Signature of the authorized person: _____

[illegible]



HERMLE Labortechnik GmbH
Siemensstraße 25
78564 Wehingen
Tel: 0 74 26-96 22-17
Fax: 0 74 26-96 22-49
Email: vertrieb@hermleLT.de
Internet: <http://www.hermle-labortechnik.de>

Technical modification rights reserved.
©HERMLE Labortechnik GmbH 2018